



INSTALLATION, OPERATION & MAINTENANCE INSTRUCTIONS

1455 Kleppe Lane ♦ Sparks, NV 89431-6467 ♦ (775) 359-4712 ♦ Fax (775) 359-7424
E-mail: haws@hawsco.com ♦ website: www.hawsco.com

No. 0002080152 (23)

Model 8317 Combination Shower Eye/Face Wash

NOTE TO INSTALLER: Please leave this information with the Maintenance Department.

LIMITED WARRANTY

HAWS warrants that this specific product is guaranteed against defective material or poor workmanship for a period of **one year from date of shipment**. HAWS liability under this warranty shall be discharged by furnishing without charge F.O.B. HAWS Factory any goods, or part thereof, which shall appear to the Company upon inspection to be of defective material or not of first class workmanship, provided that claim is made in writing to Haws within a reasonable period after receipt of the product. Where claims for defects are made, the defective part or parts shall be delivered to the Company, prepaid, for inspection. HAWS will not be liable for the cost of repairs, alterations or replacements, or for any expense connected therewith made by the owner or his agents, except upon written authority from HAWS, Sparks, Nevada. HAWS will not be liable for any damages caused by defective materials or poor workmanship, except for replacements, as provided above. Buyer agrees that Haws has made no other warranties either expressed or implied in addition to those above stated, except that of title with respect to any of the products or equipment sold hereunder and that HAWS shall not be liable for general, special, or consequential damages claimed to arise under the contract of sale.

The emergency equipment manufactured by HAWS is warranted to function if installation and maintenance instructions provided are adhered to. The units also must be used for the purpose for which they were intended. This product is intended to supplement first-aid treatment. Due to widely varying conditions, Haws cannot guarantee that the use of this emergency equipment will prevent serious injury or the aggravation of existing or prior injuries.

NO OTHER WARRANTIES EXPRESSED OR IMPLIED ARE AUTHORIZED, PROVIDED OR GIVEN BY HAWS.

SHOULD YOU EXPERIENCE DIFFICULTY WITH THE INSTALLATION OF THIS MODEL PLEASE CALL:

TECHNICAL SUPPORT: 1-800-766-5612

FOR CUSTOMER SERVICE: 1-888-640-4297

LOCATION OF UNIT: The Model 8317 Emergency Shower should be installed in close proximity to potential accident areas. It should be clearly identified, free from obstructions and easy to access.

SUPPLY LINE: The minimum recommended line size is 1-1/4" IPS with 30-90 psi (2-6 ATM) flowing pressure. Where sediment or mineral content is a problem, an inlet filter is recommended.

PLUMBING CONNECTIONS: Inlet supply is female 1-1/4" IPS.


INSTALLATION PROCEDURE

Refer to Installation Drawing for clarification. Check parts list to ensure all required parts are included. Use pipe joint sealant PST #565 (Loctite) on all connections.

- STEP 1:** Prepare the emergency shower foundation by using the floor flange assembly as a template. Locate the three holes for the 3/8-16 UNC x 6" anchor bolts (not supplied). They should extend 1" above the floor level.
- STEP 2:** Once the concrete foundation is ready (dry), assemble the floor flange assembly to the anchor bolts, then securely fasten with 3/8-16 UNC Hex Nuts and 3/8" washers (not supplied). Seal and tighten floor flange and (1-1/4") coupler. Make sure foot control holes are facing side to side.
- STEP 3:** Assemble plugged nipple (1-1/4" x 24") to the floor flange assembly and tighten with a pipe wrench. (Plugged end up).
- STEP 4:** Assemble the eyewash assembly and tee (1-1/4" x 1-1/4" x 1/2") to the 24" nipple from Step 3.
- STEP 5:** Attach nipple (1-1/4" x 31") and tee (1-1/4") to assembly and tighten.
- STEP 6:** Attach nipple (1-1/4" x 24") and tee (1-1/4" x 1-1/4" x 1) and tighten.
- STEP 7:** Attach valve assembly with nipple (1" x 9") to tee (1-1/4" x 1-1/4" x 1) with a spud wrench.
- STEP 8:** Assemble the street elbow (1") onto nipple (1" x 12-1/2").
Attach nipple onto valve.
Attach the showerhead onto the street elbow (1").
- STEP 9:** Assemble the sign onto the unit. Use the strap, screws, washers, and nuts provided, to mount the sign onto the upright. The sign should be positioned for maximum visibility.
- STEP 10:** Insert the galvanized plug into the optional supply line tee that is not used. Before connecting the supply water line to the emergency shower, **FLUSH THE LINE THOROUGHLY.**

TROUBLESHOOTING	
PROBLEM	REPAIR CHECKLIST
1. No flow.	1. Check the main shut-off valve.
2. Insufficient water flow, shower and eyewash.	2. Verify minimum 30 psi (2 ATM) flowing supply line pressure.
3. Insufficient water flow on eye/face wash, however, shower is sufficient.	3. a. Verify minimum 30 psi (2 ATM) flowing supply line pressure. b. Probable clogging of flow control due to inadequate line flushing. Unscrew eye/face wash head and remove the four screws to disassemble the head. Clean flow control and reassemble head. See SP65 Installation Drawing for details.
4. Eye/face wash stream does not meet desired eye level.	4. a. Possible blocked flow control, see above solution. b. Possible non-leveled eye/face wash assembly.
5. Insufficient water flow at the showerhead, however, eye/face wash is sufficient.	5. Probable clogging of flow control inside shower head. Unscrew shower head off of elbow and clean rubber flow control.
6. Water does not drain properly.	6. Check your drain system for debris. Check the main waste line of your building to see if it does handle the capacity required for the entire drainage system.
7. Water leak between joints or fittings.	7. Tighten loose connections. If leak persists, the application of hot water will sometimes reseal improperly applied joint compound. After 30 minutes wait and retest, a persistent leak will necessitate reassembly and additional sealant.

For more information on Haws Products, please see our website: www.hawasco.com

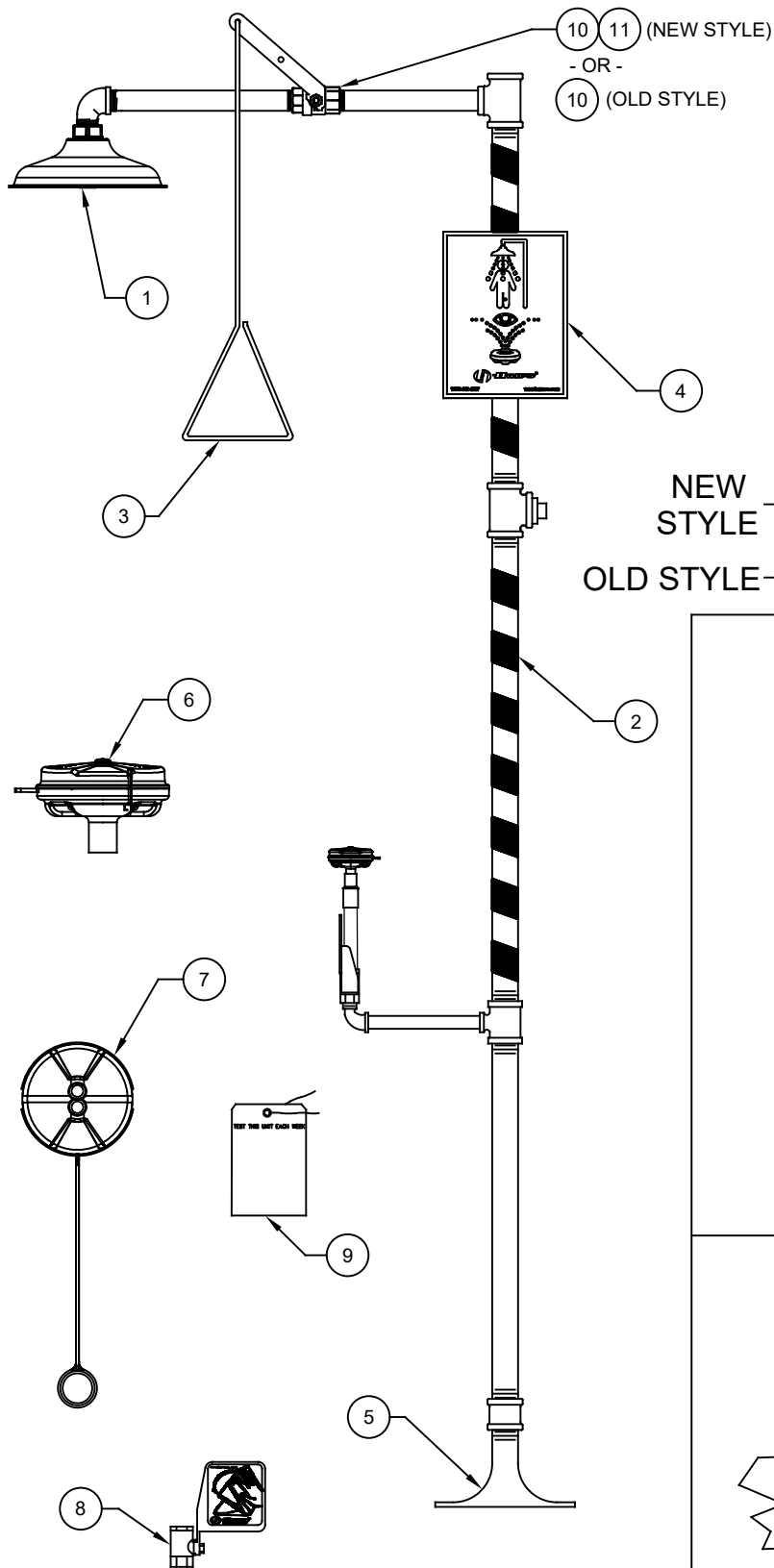


EYEWASH VALVE
ASSEMBLY

PARTS BREAKDOWN


WHEN ORDERING PARTS PLEASE SPECIFY MODEL NUMBER.

ITEM	DESCRIPTION	PART NO.
1	SHOWER HEAD	SP829
2	STRIPE TAPE	SP185
3	PULL ROD	SP200
4	SHOWER SIGN	SP178
5	FLOOR FLANGE	SP75
6	HEAD ASSY SP65	0002953160
7	DUST COVER F/SP65	9091
8	EYEWASH VALVE ASSM	SP238
9	TEST TAG	SP170

10	VALVE, 1" BALL	SP259
11	PULL ROD HANDLE	0410000858
10	VALVE/HANDLE ASSM	SP260

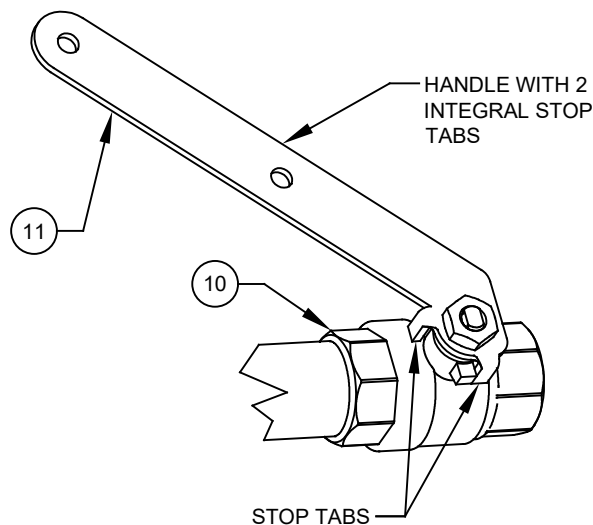

NEW STYLE VALVE & HANDLE

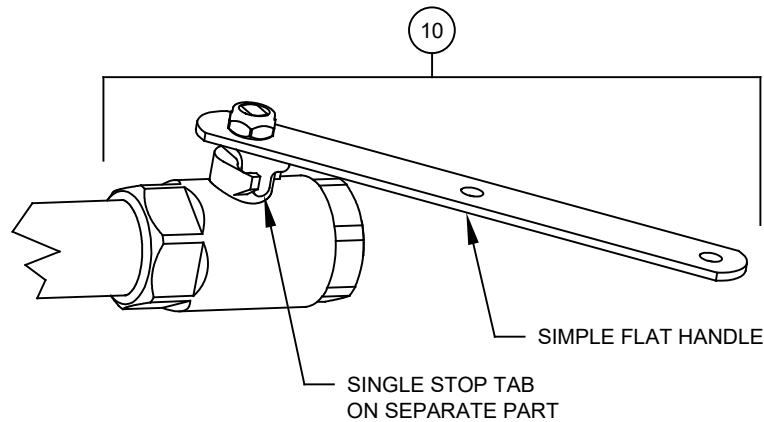


OLD STYLE VALVE & HANDLE




Haws®

1455 KLEPPE LANE
SPARKS, NEVADA 89431
(775) 359-4712 FAX (775) 359-7424
E-MAIL: HAWS@HAWS.CO.COM
WEBSITE: WWW.HAWS.CO.COM


ECN NO.	REVISED PER	BY	MODEL(S)	PART NUMBER
DRAWN	CLF	DATE	CHKD	BJM
APPROVED	JL	DATE	01/12/23	SCALE: 1:1
8317	SHOWER EYE/FACE WASH COMBINATION	REVISION	23	1

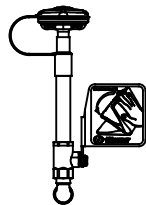

EYE/FACE WASH DETAIL

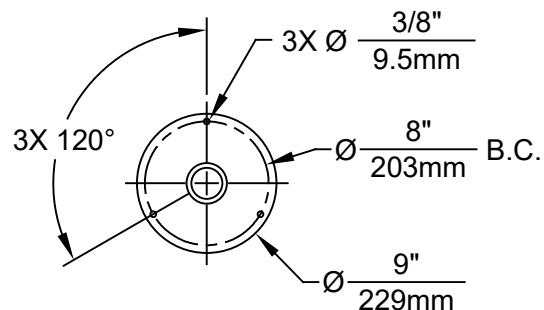

MOUNTING DETAIL

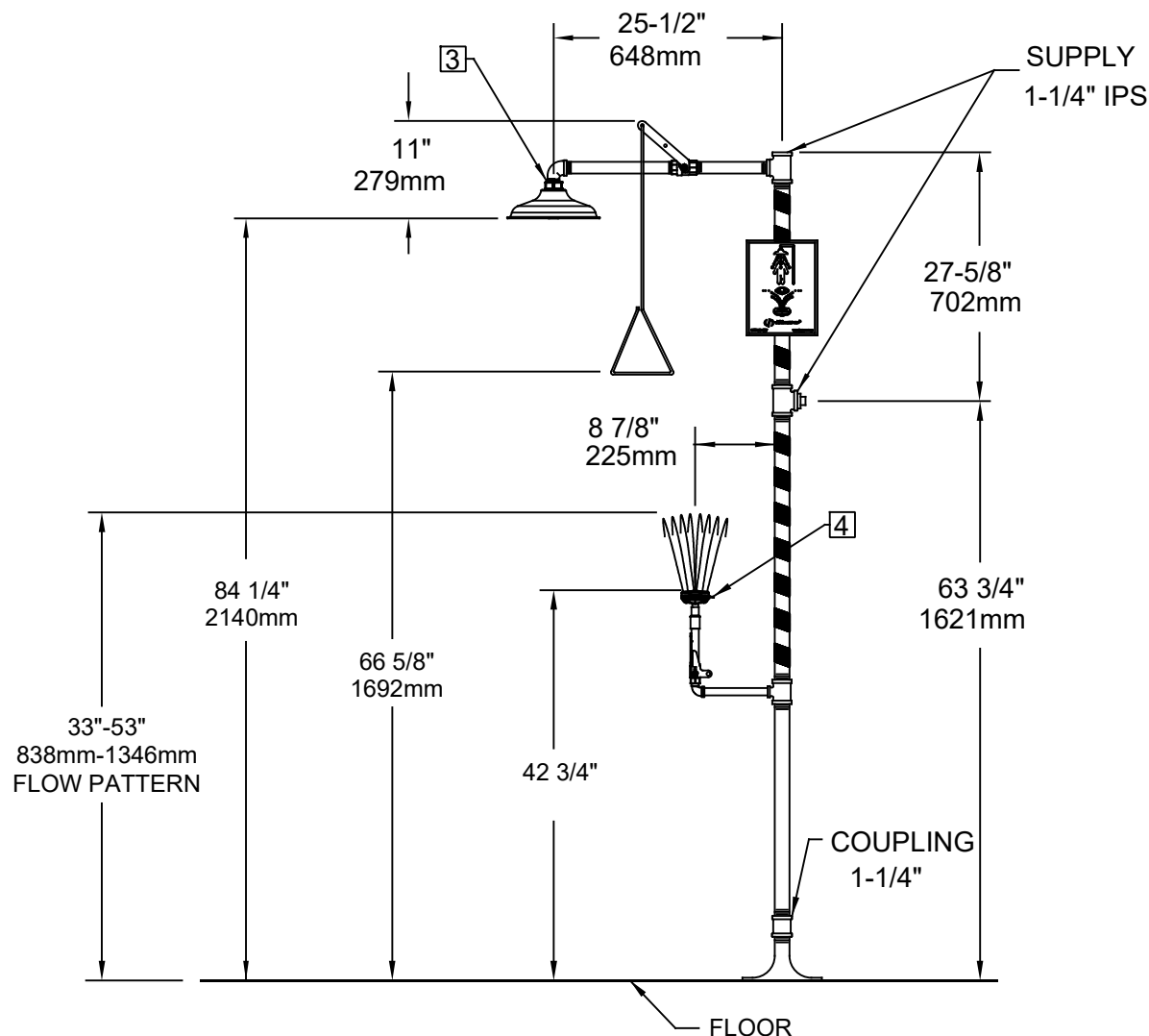

NOTES:

1. WHEN INSTALLING THIS UNIT, LOCAL, STATE, OR FEDERAL CODES SHOULD BE ADHERED TO FOR INSTALLATION DIMENSIONS (LOCATIONS), WASTE AND SUPPLY REQUIREMENTS OTHER THAN SHOWN.
2. TO COMPLY WITH ANSI Z358.1-2014 FOR EMERGENCY EYEWASH OR EYE/FACE WASH:
 - EYEWASH OR EYE/FACE WASH SHALL BE POSITIONED SUCH THAT FLUSHING FLUID PATTERN IS NOT LESS THAN 33" (838mm), NO GREATER THAN 53" (1346mm) FROM SURFACE ON WHICH USER STANDS, AND 6.0" (153mm) MINIMUM FROM WALL OR NEAREST OBSTRUCTION.
 - SHOWER SHALL BE POSITIONED SUCH THAT SHOWER HEAD IS NOT LESS THAN 82" (2083mm) AND NO GREATER THAN 96" (2438mm) FROM SURFACE ON WHICH USER STANDS.
 - SHOWER SPRAY PATTERN SHALL HAVE A MINIMUM DIAMETER OF 20" (508mm) AT 60" (1524mm) ABOVE SURFACE ON WHICH USER STANDS. CENTER OF SPRAY PATTERN SHALL BE LOCATED AT LEAST 16" (406mm) FROM ANY OBSTRUCTION.
3. DO NOT USE REGULAR PIPE SEALANT. IF NECESSARY USE DOW CORNING 832 SEALANT PLACING 5 SMALL BEADS ON MALE THREADS, JUST ABOVE THE FIRST THREAD.
4. ORIENT TAB ON EYE/FACE WASH TOWARD BACK OF SHOWER.
5. DIMENSIONS MAY VARY BY $\pm 1/2$ " (13mm).

		1455 KLEPPE LANE SPARKS, NEVADA 89431 (775) 359-4712 FAX (775) 359-7424 E-MAIL: HAWS@HAWSO.COM WEBSITE: WWW.HAWSO.COM	
ECN NO.	REVISED PER	BY:	MODEL(S)
CH	DATE	CHK'D:	8317
05/91	01/12/23	BJM	SHOWER EYE/FACE WASH COMBINATION
APPROVED:	DATE:	SCALE:	DRAWING TYPE:
JL	01/12/23	1:1	INSTALLATION
PART NUMBER		REVISION	
0002080152.D		23	
SIZE: A		SHEET 1 OF 1	