

## NOTES:

- 1. WHEN INSTALLING THIS UNIT, LOCAL, STATE, OR FEDERAL CODES SHOULD BE ADHERED TO FOR INSTALLATION DIMENSIONS (LOCATIONS), WASTE AND SUPPLY REQUIREMENTS OTHER THAN SHOWN.
- TO COMPLY WITH ANSI Z358.1-2014 FOR EMERGENCY EYEWASH OR EYE/FACE WASH:
  EYEWASH OR EYE/FACE WASH SHALL BE POSITIONED SUCH THAT FLUSHING FLUID PATTERN IS NOT
  LESS THAN 33" (838mm) AND NO GREATER THAN 53" (1346 mm) FROM SURFACE ON WHICH USER
  STANDS; AND 6.0" (153mm) MINIMUM FROM WALL OR NEAREST OBSTRUCTION.
- 3. UNIT MUST BE LEVEL FROM FRONT TO BACK AND SIDE TO SIDE
- 4 ALL THREE MOUNTING HOLES ON MOUNTING BRACKET MUST BE USED TO PROPERLY SECURE EYEWASH UNIT TO WALL.
- 5. \*\*UNIT MUST BE FILLED WITH WATER BEFORE CONNECTING TO POWER SOURCE.
- 6. PRIOR TO SERVICING/CLEANING: UNPLUG POWER CORD ON UNIT FROM POWER SOURCE AND REMOVE UNIT FROM MOUNTING BRACKET.
- 7. \*DIMENSIONS MAY VARY ±1/2" (12.7mm)

## PORTABLE EYEWASH WITH HEATER

